



Illumina Proactive Service connectivity guide

Easily connect your MiniSeq™ Sequencing System

To reduce unplanned downtime, minimize run failures and lost samples, plus resolve issues faster if and when they arise, simply follow the steps below.

Follow these steps to get connected

- From the Welcome Screen, select **Manage instrument**
- From the Manage Instrument Screen, select **System configuration**
- From the System Configuration Screen, select **Analysis configuration**
- Select **Send instrument performance data to Illumina** from the **BaseSpace™** tab
- Click **Save**

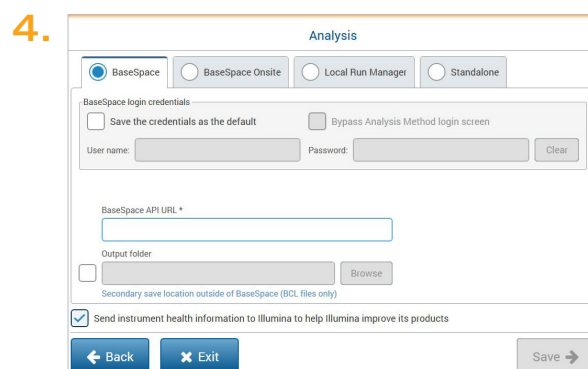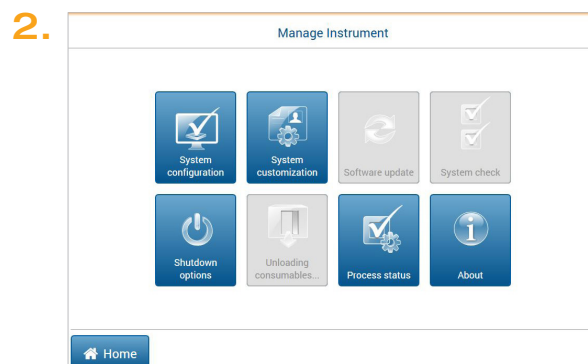

For more information, contact your local field applications scientist or field service engineer.

www.illumina.com

For Research Use Only. Not for use in diagnostic procedures.